



Table of Contents

(Use CTRL + click to follow link)

<u>Analysis Category</u>	<u>Page</u>
1. Organic Chemistry Containers	2
2. Inorganic Chemistry Containers	10
a. Radiation Containers	10
b. General Chemistry Containers	12
c. Microbiology Containers	18
d. Metals Container	19

Organic Sample Containers:




Description: 40mL Clear Glass Vial; VOC Chlorinated Drinking Water

Cap: Gray, Teflon - lined, with Septa hole

Preservation: Pre-preserved with 25 mg of Ascorbic Acid; Preserved at collection with 3 drops of Hydrochloric Acid from the blue capped, plastic dropper provided

ID Sticker: Orange dot with Ascorbic Acid label, White MDH label with analysis, expiration date and preservative information

Reference Method(s): EPA 524.2

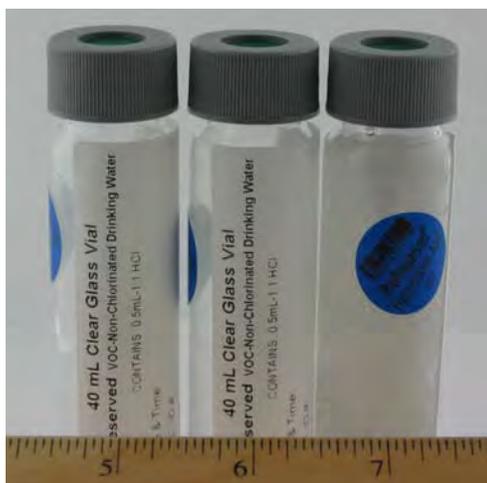

Description: 40mL Clear Glass Vial; VOC Non-Chlorinated Drinking Water

Cap: Gray, Teflon - lined, with Septa hole

Preservation: Pre-preserved with 0.5mL of Hydrochloric Acid

ID Sticker: Blue dot with HCl preservative label, White MDH label with analysis, expiration date and preservative information

Reference Method(s): EPA 524.2




Description: 40mL Clear Glass Vial; VOC Non-Potable Water

Cap: White, Teflon - lined, with Septa hole

Preservation: Pre-preserved with 0.5mL of Hydrochloric Acid

ID Sticker: Blue dot with HCl preservative label, White MDH label with analysis and preservative information

Reference Method(s): SW 8260, WI DNR, EPA 8260B

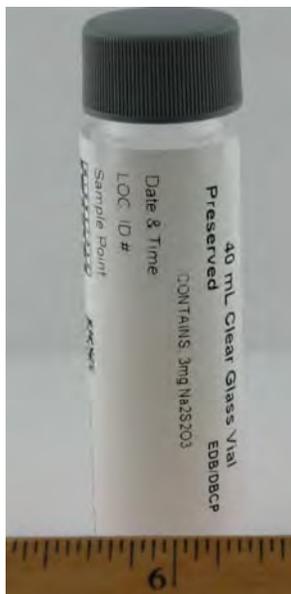


Description: 40mL Clear Glass Vial; EDB/DBCP

Cap: Gray, Teflon - lined

Preservation: Pre-preserved with 3mg of Sodium Thiosulfate

ID Sticker: White MDH label with analysis, expiration date and preservative information

Reference Method(s): EPA 504.1

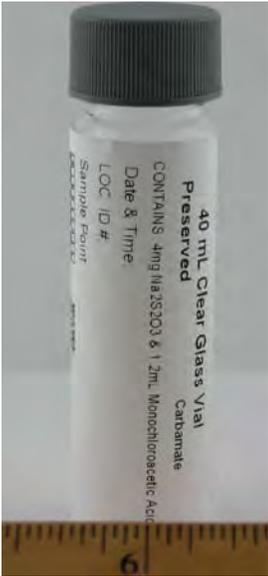

Description: 40mL Clear Glass Vial; Carbamate

Cap: Gray, Teflon - lined

Preservation: Pre-preserved with 1.2mL of Monochloro Acetic Acid and 3.2 mg of Sodium Thiosulfate

ID Sticker: Orange dot, White MDH label with analysis, expiration date and preservative information

Reference Method(s): EPA 531.1




Description: 40mL Amber Glass Vial; Glyphosate

Cap: Gray, Teflon - lined

Preservation: Pre-preserved with 4 mg of Sodium Thiosulfate

ID Sticker: White MDH label with analysis, expiration date and preservative information

Reference Method(s): EPA 547

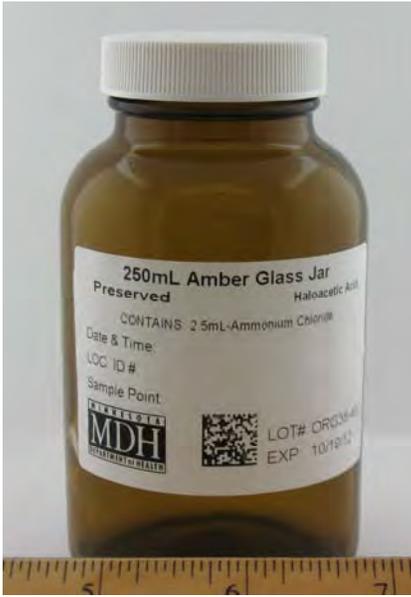


Description: 250mL Amber Glass Jar; Haloacetic Acid

Cap: White, Teflon - lined

Preservation: Pre-preserved with 2.5 mL of Ammonium Chloride

ID Sticker: White MDH label with analysis, expiration date and preservative information

Reference Method(s): EPA 552.2

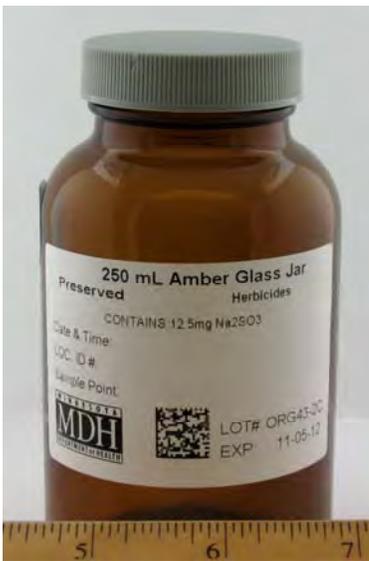


Description: 250mL Amber Glass Jar; Herbicides

Cap: Gray, Teflon - lined

Preservation: Pre-preserved with 12.5 mg of Sodium Sulfite

ID Sticker: Blue dot, White MDH label with analysis, expiration date and preservative information

Reference Method(s): EPA 515.4

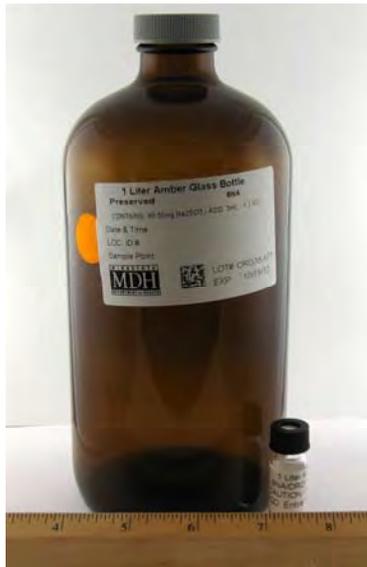

Description: 1L Amber Glass Bottle; BNA

Cap: Gray, Teflon - lined

Preservation: Pre-preserved with 40-50mg of Sodium Sulfite; Preserved at collection with 3mL of Hydrochloric Acid; add entire contents of black capped, glass vial provided

ID Sticker: Orange dot, White MDH label with analysis, expiration date and preservative information

Reference Method(s): EPA 525.2/508.1



Description: 1L Amber Glass Bottle; SVOC, PCB

Cap: Gray, Teflon - lined

Preservation: None

ID Sticker: White MDH label with analysis, expiration date and preservative information

Reference Method(s): SW 8270 D, SW 8082



Description: 1L Amber Glass Bottle; DRO

Cap: Gray, Teflon - lined

Preservation: Preserved at collection with 3mL of Hydrochloric Acid; add entire contents of black capped, glass vial provided

ID Sticker: White MDH label with analysis, expiration date and preservative information

Reference Method(s): WI DNR



Description: 60mL Clear Glass Jar; Soil VOC, GRO

Cap: White blue or gray, Teflon - lined

Preservation: Preserved at collection with 25mL of Methanol, added to jar labeled “ADD methanol,” jar labeled “NO methanol” is not preserved; add entire contents of gray capped, glass vial provided

ID Sticker: White MDH labels with fill instructions, expiration date and preservation instructions

Reference Method(s): SW 8260



Description: 60mL Clear Glass Jar; Soil DRO

Cap: White blue or gray, Teflon - lined

Preservation: None (Preserved with hexane upon receipt in the laboratory.)

ID Sticker: White MDH label with fill instructions and expiration date

Reference Method(s): WI DNR



Description: 250mL Amber Glass Jar; Organic Solids

Cap: White or gray, Teflon - lined

Preservation: None

ID Sticker: White MDH label with analysis, expiration date and preservative information

Reference Method(s): SW 8270 D, SW 8082



Description: 250mL Clear Glass Tall Jar, Flashpoint

Cap: White, Teflon- lined

Preservation: None

ID Sticker: White MDH label with analysis, expiration date and preservative information

Reference Method(s): ASTM D 3278-78



Description: 250mL HDPE Bottle; PFC

Cap: Clear

Preservation: None

ID Sticker: White MDH label with analysis, expiration date and preservative information

Reference Method(s): MDH

Inorganic Sample Containers:

Radiation Sample Containers:




Description: 4L HDPE Jug; Radiochemical Parameters Non-Potable Water

Cap: Clear

Preservation: None

ID Sticker: None

Reference Method(s): EPA 900.0, EPA 600/4-75-008, EPA 905.0, SM 7120

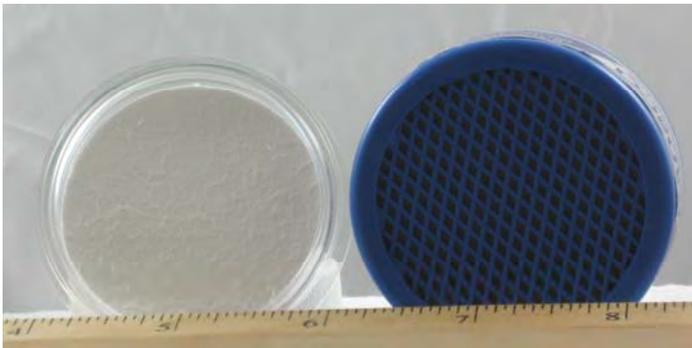

Description: 47mm Glass Fiber Filter; Radiological Parameters Air

Cap: None, filter is housed in petri dish upon sample collection

Preservation: None

ID Sticker: None

Reference Method(s): SM 7120, MDH



Description: 40mL Clear Glass Vial; Radon

Cap: Green, with septa hole

Preservation: None

ID Sticker: White MDH label with analysis, expiration date and preservative information

Reference Method(s): SM 7500-Rn



Description: 4L LDPE Cubitainer; Radiation Chemistry Parameters

Cap: White - lined

Preservation: None

ID Sticker: White MDH label with analysis, expiration date and preservative information

Reference Method(s): EPA 200.8, EPA 900.0, EPA 520/5-84-006, EPA 600/4-75-008,
EPA 905.0, EPA 903.0, EPA 904.0, SM 7120

General Chemistry Sample Containers:




Description: 1L Amber Glass Bottle; Chlorophyll-a

Cap: Green - lined

Preservation: None

ID Sticker: White MDH label with analysis, expiration date and preservative information

Reference Method(s): SM 10200 H

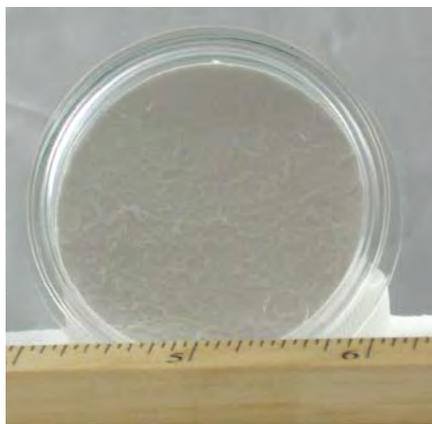


Description: 47mm Glass Fiber Filter; Chlorophyll-a, Pheophytin-a

Cap: None, filter is housed in a clear petri dish

Preservation: None

ID Sticker: None

Reference Method(s): SM 10200 H

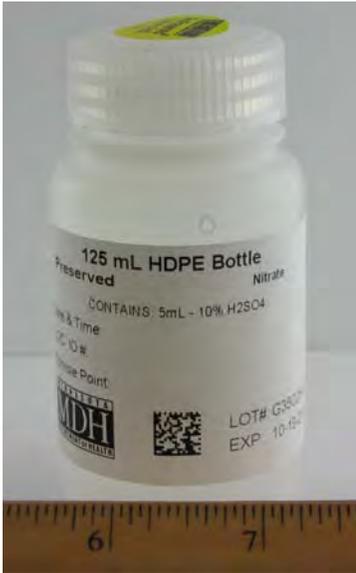

Description: 125mL HDPE Bottle; Nitrate

Cap: Clear

Preservation: Pre-preserved with 5.0mL of 10% Sulfuric Acid

ID Sticker: Yellow pre-preserved sticker, White MDH label with analysis, expiration date and preservative information

Reference Method(s): EPA 353.2, EPA 350.1




Description: 125mL HDPE Bottle; Nitrate

Cap: White - lined

Preservation: None

ID Sticker: White MDH label with analysis, expiration date and preservative information

Reference Method(s): EPA 353.2

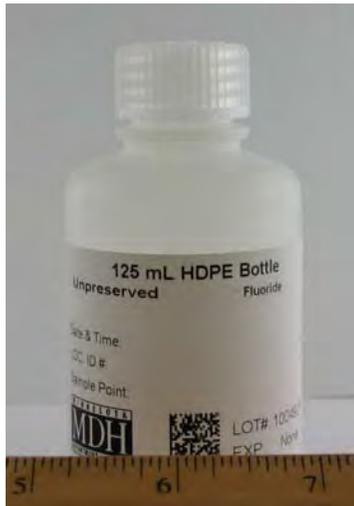

Description: 125mL HDPE Bottle; Fluoride

Cap: Clear

Preservation: None

ID Sticker: White MDH label with analysis, expiration date and preservative information

Reference Method(s): SM 4500-F C



Description: 500mL HDPE Bottle; Cyanide

Cap: White - lined

Preservation: Preserved at collection with 20mL of 2.5N Sodium Hydroxide; add entire contents of small opaque bottle provided

ID Sticker: White MDH label with analysis, expiration date and preservative information

Reference Method(s): SM 4500-CN F, EPA 335.4



Description: 250mL HDPE Bottle; Nutrient Parameters

Cap: White - lined

Preservation: Preserved at collection with 5.0mL of 10% Sulfuric Acid; add entire contents of yellow capped, plastic vial provided

ID Sticker: White MDH label with analysis, expiration date and preservative information

Reference Method(s): EPA 365.1, EPA 350.1, EPA 351.2, EPA 353.2, EPA 410.4, SM 5310 C




Description: 125mL HDPE Bottle; General Chemistry Parameters

Cap: White - lined

Preservation: None

ID Sticker: White MDH label with analysis, expiration date and preservative information

Reference Method(s): SM 4500 NO2 B, SM 2320 B, SM 4500-F C, SM 4500-SiO2 C, SM 4500 PE, EPA 365.1, EPA 300.1

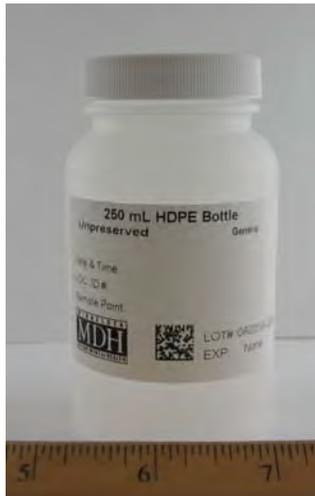


Description: 250mL HDPE Bottle; General Chemistry Parameters

Cap: White - lined

Preservation: None

ID Sticker: White MDH label with analysis, expiration date and preservative information

Reference Method(s): SM 5910 B, SM 5310 C, EPA 365.1, EPA 300.1

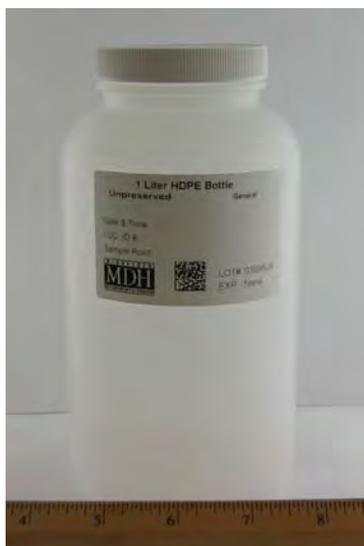

Description: 1L HDPE Bottle; General Chemistry Parameters

Cap: Clear

Preservation: None

ID Sticker: White MDH label with analysis, expiration date and preservative information

Reference Method(s): SM 2540 B, SM 2540 C, SM 2540 D, SM 2540 E, SM 2130 B,
SM 4500-H+B, SM 2510 B, SM 2320 B, SM 4500-SiO₂ C, SM
5910 B, SM 2120 B, EPA 300.1, EPA 160.4, EPA
365.1



Description: 2L HDPE Bottle; General Chemistry Parameters

Cap: Clear

Preservation: None

ID Sticker: White MDH label with analysis, expiration date and preservative information

Reference Method(s): SM 2540 B, SM 2540 C, SM 2540 D, SM 2540 E, SM 2130 B, SM 4500-H+B, SM 2510 B, SM 2320 B, SM 4500-SiO₂ C, SM 5910 B, SM 4500-P E, EPA 300.1, EPA 160.4, EPA 365.1,



Description: 250mL HDPE Jar; Inorganic Solids

Cap: White


Preservation: None single container may be used for soil analysis of both metals and chemistry parameters

ID Sticker: White MDH label with analysis, expiration date and preservative information

Reference Method(s): Quik Chem 13-115-01-1-B, Quik Chem 13-107-06-2-D, EPA 9045 D, EPA 300.1



Microbiology Sample Containers:

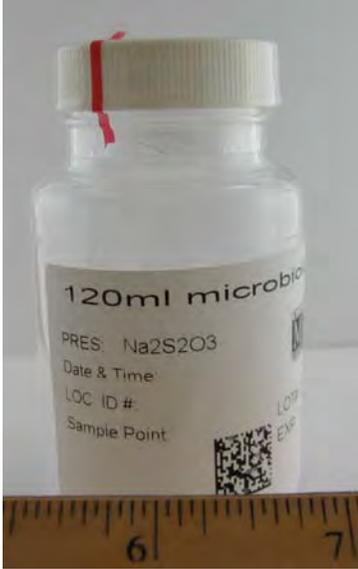

Description: 120mL HDPE (Sterile) Bottle; Microbiological Parameters

Cap: White - lined

Preservation: Pre-preserved with 10–35mg of Sodium Thiosulfate

ID Sticker: White MDH label with analysis, expiration date and preservative information

Reference Method(s): EPA 1103.1, SM 9222 B, SM 9222 G2, SM 9222 D, SM 9230 C, SM 9215 B, SM 9213 E, SM 9223

Metals Sample Containers:




Description: 250mL HDPE Jar; Inorganic Solids

Cap: White

Preservation: None; single container may be used for soil analysis of both metals and chemistry parameters

ID Sticker: White MDH label with analysis, expiration date and preservative information

Reference Method(s): SW 6010 B, SW 6020, SW 7471, MDH

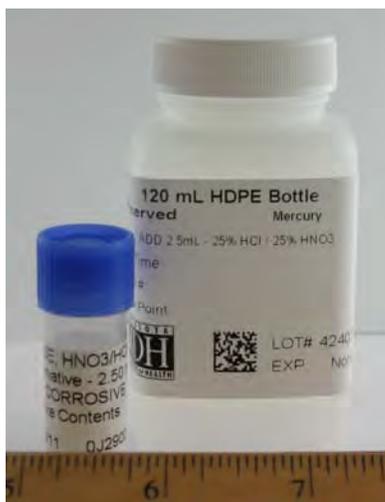

Description: 120mL HDPE Bottle; Mercury

Cap: White - lined

Preservation: Preserved at collection with 2.5 mL of 25% Hydrochloric Acid and 25% Nitric Acid; add entire contents of blue capped, plastic vial provided

ID Sticker: White MDH label with analysis, expiration date and preservative information

Reference Method(s): EPA 245.2, SW 6020



Description: 250mL HDPE Bottle; Metals Parameters

Cap: White - lined

Preservation: Preserved at collection with 2.5mL of 20% Nitric Acid; add entire contents of red capped, plastic vial provided

ID Sticker: White MDH label with analysis, expiration date and preservative information

Reference Method(s): SM 2340 B, EPA 200.7, EPA 200.8, SW 6020